

Cellular Imaging

Research Administration
Seattle, WA • 501(c)(3) Nonprofit



Fred Hutch's Shared Resources are catalysts for lifesaving discoveries. This uniquely centralized program of 15 specialized core facilities and scientific services drives advances by integrating dedicated experts and cutting-edge technologies across the entire research pipeline, from basic science to clinical trial.

Nikon TIRF/STORM

Widefield fluorescence and total internal reflection fluorescence microscope with super resolution (N-STORM)

Excitation sources

- Lasers: 405, 488, 561, 640 and 640 nm high power for STORM imaging

Objectives

- 10x/0.3, 60x/1.49, 100x/1.49

Camera

- Andor iXon X3 EMCCD camera with 512 x 512 pixel chip (16 micron pixels)

Capabilities

- Four-channel (blue, green, red, far red) TIRF
- Super-resolution localization microscopy (N-STORM)
- Time-lapse imaging

Recommended uses

- Imaging and quantitation of radioactivity, fluorescence and chemiluminescence
- Gel and blot imaging
- Alternative to fluorescence plate reader
- DNA gels (Southern blots)
- Western blots

General information

Total internal reflection fluorescence, or TIRF, microscopy uses evanescent wave technology to restrict fluorescence excitation to a very thin layer of approximately 100 to 200 nm near the coverslip surface, greatly reducing background fluorescence and improving image contrast. Therefore, TIRF is ideally suited for the detection of faint signals at or near the cell surface, and for the detection of molecules in in vitro reconstituted systems. Stochastic optical reconstruction microscopy, or STORM, is a super-resolution technique that relies on the precise localization of individual molecules to provide image resolution down to a few tens of nanometers.

Our Nikon system combines TIRF and STORM on a single microscope. The system is equipped with the latest EMCCD technology (Andor iXon X3 EMCCD camera), capable of providing single-molecule detection sensitivity. The system includes a four-laser launch with 405, 488, 561 and 640 nm laser lines for excitation of blue, green, red and far red dyes and fluorescent proteins. The system is capable of TIRF (as well as conventional fluorescence) imaging in four channels (blue, green, red, far red).

The microscope is built on a Nikon Ti fully automated inverted microscope with Perfect Focus, motorized x,y stage and fast piezo z stage. It uses Nikon NIS Elements software.

The Nikon TIRF/STORM microscope was funded in part by a generous grant by the M.J. Murdock Charitable Trust.

LEARN MORE

Cellular Imaging Core
206.667.4205
imaging@fredhutch.org

