

Glacios checklists

What to expect when you arrive

<input type="checkbox"/>	K3 computer:
<input type="checkbox"/>	Digital Micrograph
<input type="checkbox"/>	Notepad: list of grids in microscope
<input type="checkbox"/>	Photo Viewer: atlases of all grids (jpgs)
<input type="checkbox"/>	Remote Desktop: signed into WARP PC
<input type="checkbox"/>	Glacios computer
<input type="checkbox"/>	Tecnai User Interface
<input type="checkbox"/>	TIA
<input type="checkbox"/>	SEM server

Glacios startup

<input type="checkbox"/>	Start session on iLab Kiosk (support PC)
<input type="checkbox"/>	Check vacuum readings (Setup tab)
<input type="checkbox"/>	Check nitrogen levels (Autoloader tab)
<input type="checkbox"/>	Select "Turbo Auto Off" (Autoloader tab)
<input type="checkbox"/>	Insert 50um C2 aperture
<input type="checkbox"/>	Lower the flu screen
<input type="checkbox"/>	IMAGE! (go to leginon protocol)

Glacios shutdown (without collection)

<input type="checkbox"/>	Label grids to save/discard
<input type="checkbox"/>	Close column valves

Glacios shutdown (with collection)

<input type="checkbox"/>	Label grids to save/discard
<input type="checkbox"/>	Make sure SerialEM, transfer, and WARP are running
<input type="checkbox"/>	Check email at end on SerialEM
<input type="checkbox"/>	Check column valves close SerialEM
<input type="checkbox"/>	"Fill Now" (Autoloader tab) before you leave for the night